

Data Collection Software Supplemental Maintenance Release Software Patch CE_DC_300-002 – Part Number 4375632

1 Data Collection Software v3.0 – Communication Issue

Some Applied Biosystems 3130 /3130*xl* Genetic Analyzer and 3730/3730*xl* DNA Analyzer users have reported a communication problem between the Data Collection Software and Sequencing Instrument. Some of the symptoms include:

- Unable to perform a run since the Green Start button is disabled
- Unable to perform a Spatial Calibration since the “Start” Spatial Calibration button is disabled

2 Data Collection Software versions affected

This software patch is specifically designed for:

- 3130 and 3130*xl* Data Collection Software v3.0
- 3730 and 3730*xl* Data Collection Software v3.0

3 Patch Details

This software patch modifies the Data Collection Software startup process to automatically delete a set of corrupted temporary files.

4 Patch Instructions

1. Stop the Data Collection Software.
2. Copy the patch installer file to a directory on your computer.
3. Launch the patch by double-clicking on the file icon.
4. Follow the on-screen directions in the installation wizard to complete the patch installation.

5 Disclaimer

This Supplemental Maintenance Release Software Patch (SMR) has been tested for specific functionality but has not been through a complete regression test. Due to the nature of SMRs, Applied Biosystems makes no representations regarding their use or performance and accepts no liability as a result of the use of SMR software. This SMR is provided "AS IS". APPLIED BIOSYSTEMS PROVIDES NO WARRANTIES EITHER EXPRESS OR IMPLIED, INCLUDING IMPLIED WARRANTIES OF MERCHANTABILITY AND FITNESS FOR A PARTICULAR PURPOSE ALL OF WHICH ARE HEREBY DISCLAIMED. IN NO EVENT WILL APPLIED BIOSYSTEMS BE LIABLE FOR ANY INDIRECT, INCIDENTAL, CONSEQUENTIAL, SPECIAL OR EXEMPLARY DAMAGES, INCLUDING, BUT NOT LIMITED TO, LOSS OF PROFITS, OR LOSS OF BUSINESS OPPORTUNITY, IN RELATION TO THE PERFORMANCE, USE, OR OPERATION OF THIS SMR, EVEN IF SUCH DAMAGES ARE FORESEEABLE AND WHETHER OR NOT APPLIED BIOSYSTEMS HAS BEEN ADVISED OF THE POSSIBILITY THEREOF. The customer shall have sole responsibility for adequate protection and back-up of data used in conjunction with this SMR.

For Research Use Only. Not for use in diagnostic procedures.

Applied Biosystems and AB (Design) are registered trademarks and Applera is a trademark of Applera Corporation or its subsidiaries in the US and/or certain other countries. All other trademarks are the sole property of their respective owners.